Temperature Controlled Shaking

Leading Safety Standards

Superior Ease of Use

Reduced Cost of Ownership

Operating Manual

Platform Shaker Hei-MIX Titramax 100/101/1000,
Vibramax 100/101 and Rotamax 120.

Operating manual must be read before initial start-up.

Please follow the safety instructions provided.

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Seite 26

About this Document	4
Version and variants	4
About this manual	4
Terms and signs used.....	5
Safety Instructions	6
General safety instructions.....	6
EC Declaration of Conformity.....	6
Intended use.....	6
Installation / Electrical safety	7
Personnel qualification.....	7
Operating company's obligations.....	7
Installation site	7
Changes to the unit	7
Safety for the personnel	8
Safety during use	8
Disposal	8
Device Description	9
Device overview	9
Vibramax 100 overall view	9
Vibramax 110 overall view	9
Titramax 100/101/1000 overall view	10
Rotamax 120 overall view	10
Control panel	10
Start-Up	11
Setting up device	11
Connect/disconnect power cord	11
Switch unit on and off	12
Install vessels on shaker plate.....	12
Operation	14
Operation and controls	14

Cleaning and Maintenance	15
Cleaning.....	15
Maintenance	15
Troubleshooting	16
Possible errors and how to resolve	16
Assembly	17
Electrical connections.....	17
Disassembly and Storage	18
Dismantling, transportation and storage	18
Dismantling	18
Transportation and storage	18
Accessories and Spare Parts	19
Scope of delivery.....	19
Accessories.....	19
Attachments	20
Technical data	20
Performance range / admissible load	21
Service	22
Contact / Technical Service	22
Warranty	23
Confirmation of condition of unit.....	24

> Version and variants

Version

This manual describes the function, operation and maintenance of the platform shakers Hei-MIX Titramax 100/101/1000, Vibramax 100/101 and Rotamax 120.

Version	Alteration date
0.1	10.2015

Variants

The units are available in different configurations. Certain features and functions apply to specific configurations only. The varieties are documented in this manual.

> About this manual

This operating manual is a component of the device described here.

- Please read this manual carefully and obey all safety and warning notices.
- Ensure that every operator reads this manual.
- Ensure that this manual is accessible for every operator.
- Pass on the operating manual to the subsequent owner.

For the current version of this operating manual in pdf format, please go to:

<http://www.heidolph-instruments.com/Support/Operation Manuals/Shakers & Mixers/Platform Shakers/Hei-MIX Titramax 100/101/1000, Vibramax 100/101 and Rotamax 120>

In addition, please observe the regional regulations.

➤ Terms and signs used

In this manual you will find standardized terms and signs to warn you of possible dangers or give important hints. Please take special notice to these signs to avoid personal injury or damage to the appliance.

The following terms and type of signs (basic drawing elements) are used:

Symbol	Additional term / Description
Warning signs 	The yellow triangle indicates hazardous situations. It is used in combination with the following terms: DANGER: Indicates a hazardous situation which, if not avoided, will result in irreversible injury or death. WARNING: Indicates a hazardous situation which, if not avoided, could result in serious injury or material damage. CAUTION: Indicates a hazardous situation which, if not avoided, may result in a minor injury or material damage.
Prohibitory signs 	Not to do: The red circle indicates a situation that should be avoided under all circumstances to avoid injury or damage.
Mandatory signs 	To do: The blue circle indicates important information. Please obey to avoid property damage.

Other signs used

Symbol	Description
→	Handling instruction, action required
✓	Result of action
▪ ▪ ▪	List of information
a. b. c.	List of variants

➤ **General safety instructions**

EC Declaration of Conformity

This device complies with the following EC-Directives*:

- 2006/42/EC Machinery Directive
- 2006/95/EC Low-Voltage Directive
- 2004/108/EC Electromagnetic Compatibility Directive

*Also see attachment „EC Declaration of Conformity“.

The device has been constructed according to state-of-the-art technology and recognized safety regulations. However, risks may still arise during installation, operation and maintenance.

→ Please ensure the operating manual is available at all times.

The device may only be used under the following circumstances:

- Only operate the device, if it is in full working order.
- Ensure all operators of the device possess the necessary safety and risk awareness.
- Operate the device according to the instructions stipulated in this manual only.
- If there is something you do not understand, or certain information is missing, ask your manager or contact the manufacturer.
- Do not do anything on the device without authorization.
- Only use the device according to its intended use.

Intended use

The device is intended for use by trained and authorized personnel only.

The device is suitable for the following use:	<ul style="list-style-type: none"> ▪ mixing ▪ shaking ▪ separating ▪ blending of solvents
The operation for the device is suitable in one of the following locations:	<ul style="list-style-type: none"> ▪ in any research application ▪ approved for installation in gassing and conditioning cabinets

Improper use

Any use which deviates from the device’s intended use is considered to be improper. The manufacturer does not accept liability for any damages resulting from non-permitted uses. The risk is carried by the operator alone.

Installation / Electrical safety

- The device may only be connected when the mains voltage corresponds to the information on the type plate of the unit.
- The mains connection must be accessible at all times.
- Repairs may only be performed by a qualified electrician.
- Never operate the unit with a damaged power cord.
- Always turn the unit OFF and disconnect mains power before performing any maintenance or service.

Personnel qualification

- The device may only be operated by qualified persons.
- The device may only be operated by individuals who have been instructed in its proper use by qualified persons.
- The device may only be operated and maintained by persons who are of legal age.
- Other personnel may only work with the unit under continuous supervision of an experienced qualified operator.
- This manual must be read and understood by all persons working with the device.
- The personnel must have received special safety instructions in order to guarantee responsible and safe work procedures.

➤ Operating company's obligations

Installation site

- The device must be positioned in a suitable location.
- The device must be installed sufficiently stable.
- All screw connections must be securely tightened.
- Installation and operation of the device is only permitted in facilities which are fitted with the appropriate laboratory equipment (e.g. with air extraction units).
- We recommend operation in fume hood enclosures.
- The device may only be operated in enclosed spaces and under the following environmental conditions:

Ambient temperature

- 5 - 31 °C at 80 % relative humidity, no condensation
 - 32 - 40 °C decrease linearly till max. 50 % relative humidity
-

Changes to the unit

- No unauthorized changes may be made to the unit.
- No parts may be used which have not been approved by the manufacturer.
- Unauthorized changes result in the EC Declaration of Conformity losing its validity, and the appliance may no longer be operated.
- The manufacturer is not liable for any damage, danger or injuries that result from unauthorized changes or from operating the unit other than described in this manual.

Safety for the personnel

- Ensure that only qualified personnel operate the device.
- Observe the following regulations:
 - Laboratory guidelines
 - Accident prevention regulation
 - Ordinance on Hazardous Substances
 - Other generally accepted rules of safety engineering and occupational health
 - Local regulations

➤ Safety during use

- Wear the appropriate protective clothing when working on the device (clothing, protective glasses and if necessary, safety gloves).
- Do not use the device in potentially explosive areas. The device is not protected against explosion. There is no explosion or ATEX protection available.
- Do not operate or assemble devices in the vicinity which are emission or radiation sources (electromagnetic waves) for the frequency range (3×10^{11} Hz to 3×10^{15} Hz).
- Eliminate errors immediately.
- Do not use abrasive material to clean the glass surfaces. Only wipe with damp cloths.
- Always switch the device off after use.

➤ Disposal

- Check the device components for hazardous substances and solvents.
- Clean all components before disposal.
- Dispose of the device according to the valid national legal regulations.
- Dispose of the packaging material in accordance with the valid national legal regulations.

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Analog knob: to set speed

Analog knob: to set time

Select mode switch: switch from timer to continuous mode

ON/OFF switch

> Setting up device

Warning:
Risk of slipping device!

As a result of vibrations generated during operation the device might slide off the table top.

The surface must be smooth and clean and should be properly leveled.

- Locate the shaker on a stable, horizontal surface.
- Clean surface and feet with a damp cloth and ethanol regularly.

> Connect/disconnect power cord

Connect power cord

The power cord comes with a three-pole plug and a recess on the bottom side.


Unit plug

Power cord

The socket is located on the back side of the unit.

- Connect plug to the socket with the recess facing down.
- Push plug up against socket until securely attached.
- Connect plug to power socket.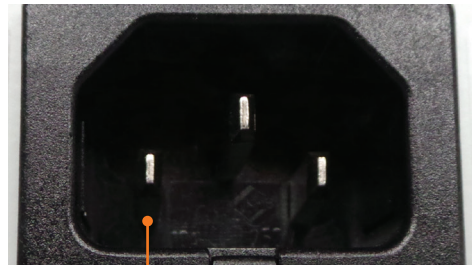

Unit socket

Disconnect power cord


- Disconnect plug from power socket first.
- Then disconnect plug from socket on unit.

➤ Switch unit on and off

Switch unit on


- Press the On/Off switch at the lower side.
- ✓ The green control light switches on.
- ✓ The unit is switched on.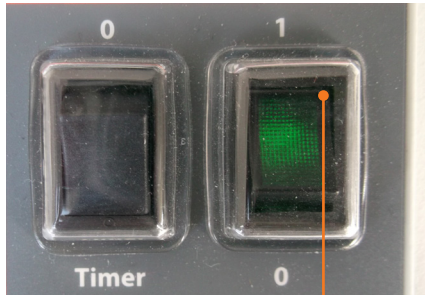

Switch On/Off: device switched on

Switch unit off

- Press the On/Off switch at the upper side.
- ✓ The green control light switches off.
- ✓ The unit is switched off.

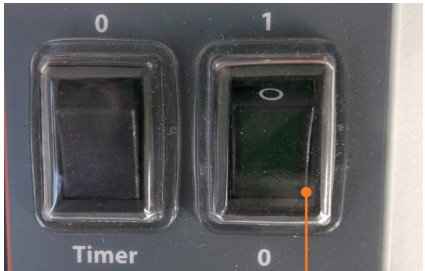Switch On/Off: device switched off

➤ Install vessels on shaker plate

Use optional attachment to secure vessels on shaker plate. Attachments with clamping rollers for a variety of flasks, beakers, and test tubes are available. Microtiter plates nest on the shaker plate, no optional holders are required.

In special cases, and at low shaking speed, e.g. Petridishes may be placed on the anti-skid rubber plate that comes with your item as a standard option.

- Before starting shaking, ensure vessels and the like are properly secured on the shaker plate.

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Microtiter plates are loaded in indentations in the rubber mat, shipped with your shaker as a standard option. They are positively locked in these indentations, the rubber mat itself can be replaced at ease.

Start-Up

VIBRAMAX 100 / ROTAMAX 120 / attachment with clamping rollers

Vessels are located on an anti-skid rubber mat, shipped with the machine. The rubber mat itself can be replaced at ease.


Vessels can be secured by an optional attachment with clamping rollers either. Standard kit includes 2 ea. clamping rollers. Extra clamping rollers can be ordered.

This attachment is installed on the shaker plate and secured with screws from your hardware bag.

VIBRAMAX 110 / attachment for test tubes

Test tube are seated on the rubber mat by hand.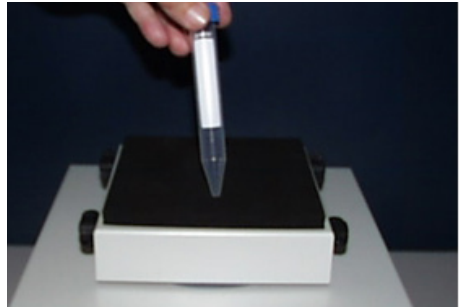

Several test tubes may be shaken at the same time, using the optional test tube attachments. These attachments are secured on the shaker table with wing bolts. Test tubes are plugged into "collets".

➤ Operation and controls

Caution:
Risk of breakage and spillage

If the vessels are not attached securely they may fall out and spill the sample.

Prior to switching the unit on ensure that all vessels are securely fastened.

Warning:
Risk of poisoning

Open vessels and too high vibration frequencies may result in samples splashing.

Close all vessels tight especially if sample is dangerous or toxic.

Adjust vibration frequency step by step until you have reached the required rpm settings..

Wear safety glasses and especially with dangerous and toxic samples appropriate protective clothing.

A single flask should be arranged in the middle of the shaker plate.

Several flasks should be distributed equally on the plate.

- ➔ Turn item ON with the On/Off switch.
- ➔ Set shaker frequency with speed setting knob.
- ➔ Select continuous mode (symbol **O**) or timer controlled mode (timer) with select mode switch.
- ✓ In the timer mode, shaking action can be selected between 0 and 120 minutes.
- ✓ After time elapsed, item will stop shaking. A buzzer sounds.
- ✓ In case of power supply failure the timer will continue running.

At high loads (load bearing capacity of shaker plate) and high frequency of shaking, always make reference to load graph as applicable for your item (refer to „Technical data“).

When using shakers inside conditioning cabinets, make reference to ambient conditions as stipulated in the „Technical data“ section.

> Cleaning

All surfaces of the unit allow for cleaning with a damp cloth and if required a mild soap lotion.

Attention:
Damage to the surfaces

All surfaces can be damaged by improper cleaning.

In any case do not use:

- Chlorine bleach or any cleaning agent including chlorine
- Ammonia
- Abrasive cleaning agents such as cleaning rags, scrubbing agents or any other agents which include metal components

As a result of vibrations generated during operation the device might slide off the table top.

To ensure a stable position clean surface and feet with a damp cloth and ethanol regularly.

> Maintenance

The unit is maintenance-free. Any necessary repair must be performed by an authorized Heidolph distributor.

Please contact Heidolph Instruments or your local Heidolph Instruments distributor.

Troubleshooting

➤ Possible errors and how to resolve

Malfunctions	Possible reason	Troubleshooting
On-/Off switch does not enlighten	No power	<ul style="list-style-type: none"> ▪ Check size and shape of plug and compatibility to your electrical socket ▪ Check circuit breakers
	Light diode defective	Contact your local Heidolph Instruments distributor
Unit stops shaking	Pre-set timer elapsed	Restart timer, if necessary
	Thermal motor circuit breaker triggered by motor overload.	Wait 20 minutes, decrease overall load
	<ul style="list-style-type: none"> ▪ Mechanical parts broken ▪ Electronic failure 	Contact your local Heidolph Instruments distributor

If you experience a malfunction which can not be resolved, please contact your authorized Heidolph Instruments distributor immediately.

> Electrical connections

- This unit must **only** be connected **by a professional electrician**.
- **Any repair** on this unit must **only** be carried out **by a qualified specialist**.

Warning: Electric shock!

This unit must only be connected to a grounded electrical socket.

Repairs are allowed by qualified professionals only. Any improper repair can result in a dangerous situation. Contact your local Heidolph Instruments distributor for any repair you may have.

The unit and voltage must match. The type label on the back side of the unit provides all voltage specifics.

When shipped, the unit is grounded. In case the original plug is changed, the new plug must have a protective conductor!

Color code for electrical connections:

Green/Yellow	PE: Protective conductor (Earth)
Blue	N: Neutral conductor
Brown	P: Phase
Black	P: Phase
Grey	P: Phase

The unit is connected with the power cord supplied to the electrical socket. It plugs into the housing on the rear side of the unit.

For countries utilizing a plug other than the standard one supplied:

- The plug supplied may only be changed by a professional electrician.
- If utilizing an adaptor ensure that it is approved by the local regulations.

➤ Dismantling, transportation and storage

Dismantling

Warning: Risk of injury!

While the unit is plugged in you might be at harm if the device is accidentally switched to ON mode.

Prior to dismantling always disconnect the plug from the mains.

- ✓ The unit is switched off.
- Disconnect plug from power socket.
- Disconnect plug from socket on the unit.

Transportation and storage

- Store and transport the unit and its components only if they were emptied and cleaned in the original packing material.
- Alternatively store and transport the unit in a container which prevents damages during transportation.
- Seal the packing carefully against unauthorized or mistaken opening.
- Store the unit in a dry place.

Improper transportation may result in damages to the system and the mechanical components.

Avoid any kind of shocks during transportation.

➤ Scope of delivery

Component	Variant	Quantity	Product number
Hei-MIX	Titramax 100	1	544-11200-00
	Titramax 101	1	544-11300-00
	Titramax 1000	1	544-12200-00
	Vibramax 100	1	544-21200-00
	Vibramax 110	1	544-31200-00
	Rotamax 120	1	544-41200-04
Operating Manual English / German		1	01-005-002-34
Warranty registration / Confirmation of condition		1	01-006-002-78
EC Declaration of Conformity		1	01-001-025-06
Power cord		1	

➤ Accessories

Component		Quantity	Product number
Vibramax 100	Attachment with 2 clamping rollers	1	549-81000-00
Rotamax 120	Clamping roller (additional)	1	11-008-007-08
Vibramax 110	Attachment for 49 test tubes, 12 mm Ø	1	549-82000-00
	Attachment for 36 test tubes, 16 mm Ø	1	549-83000-00

Additional accessories are shown in the general catalog or on our website at www.heidolph-instruments.com

➤ Technical data


Standard supply voltage	1 x 230 V (50/60 Hz)				
Operational voltage	100-240 V (50/60 Hz)				
Protection class (DIN EN 60529)	IP30				
Drive	EC-motor				
Overheat protection	self-resetting				
Speed control	analog				
Operation mode	timer / continuous				
Permissible ambient conditions	<ul style="list-style-type: none"> ▪ 5 - 40 °C ▪ 80 % relative humidity, no condensation at 5 - 31 °C, linear decreasing to 50 % at 40 °C 				
Modell Hei-MIX	Titramax 100/101	Titramax 1000	Vibramax 100	Vibramax 110	Rotamax 120
Dimensions (wxhxd) (mm)	245x125x310	320x125x375	245x125x310	245x146x310	245x125x310
Load capacity (kg)	2	5	2	2	2
Motion	circular vibration			circular vibration	rotation
Orbit (mm)	1,5 / 3	1,5	3	1,5	20
Rotation speed (rpm)	150 - 1.350			150 - 2.500	20 - 300
Supply power (W)	31			46	33
Weight (g)	5,5	6,5	5,5	5,0	5,5


Performance range / admissible load

Hei-MIX Titramax 100/101/1000


Hei-MIX Vibramax 100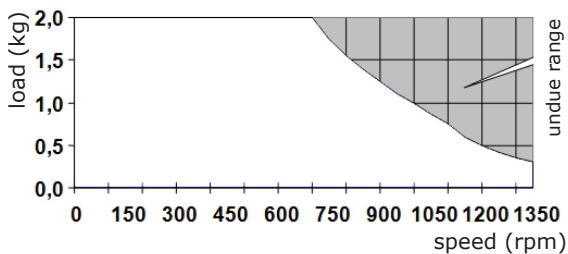

Hei-MIX Vibramax 110

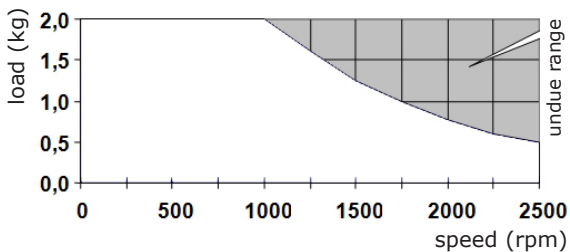Hei-MIX Rotamax 120

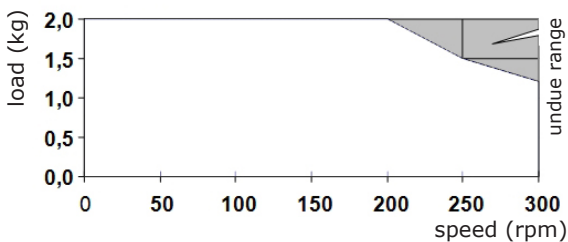➤ Contact / Technical Service

Questions / Repair work

If any aspect of installation, operation or maintenance remains unanswered in the present manual, please get in touch with the following address.

For repairs please call Heidolph Instruments or your local authorized Heidolph Instruments distributor.

Warning: Danger of poisoning!

Contaminated units can lead to severe injury or death of our employees!

When shipping items for repair that may have been contaminated by hazardous substances, please

- advise exact substance
- take adequate protective action towards our parts receiving and service personnel
- mark the pack in accordance with Ordinance on Hazardous Substances

Included at the end of this manual you will find a „Confirmation of condition of unit“.

→ Kindly copy and fill in this form and submit it prior to shipping the unit for repair.

Contact details:

Heidolph Instruments Germany

Heidolph Instruments GmbH & Co. KG
Technical Service
Walpersdorfer Str. 12
D-91126 Schwabach / Germany

Tel.: +49 – 9122 - 9920-74
Fax: +49 – 9122 - 9920-84

E-Mail: service@heidolph.de
www.heidolph.de

Heidolph Instruments North America

Phone: 1-866-650-9604
E-mail: service@heidolph.com
www.heidolphNA.com

Heidolph Instruments United Kingdom

Phone: 01799 - 5133-20
E-mail: service@radleys.co.uk
www.heidolph-instruments.co.uk

All other countries

You will find contact details of your local Heidolph distributor at www.heidolph.com

➤ Warranty

Heidolph Instruments provides a 3-year warranty for the products described here (excluding glass and wearing parts) if you register using the warranty card enclosed or online (www.heidolph.com). The warranty is valid from the point of registration. The serial number is also valid without registering.

The warranty covers part and manufacturing defects.

In the case of a part or manufacturing defect, the device shall be repaired or replaced free of charge under the terms of the warranty.

Heidolph Instruments shall not assume liability for any damages incurred as a result of improper handling or transport.

Warranty claim?

- ➔ Please inform Heidolph Instruments or your local Heidolph distributor should you wish to make a warranty claim.

➤ Confirmation of condition of unit

→ In the case of repair, copy and complete the Confirmation of condition of unit and send it to Heidolph Instruments.

1. Details about the unit

Product number _____

Serial number _____

Reason for repair _____

2. Has the device been cleaned, decontaminated/sterilized?

Yes _____ No _____

3. Is the unit in a condition which does not represent any health threats for the staff of our service department?

Yes _____ No _____

If not, which substances has the unit come into contact with?

4. Legally binding declaration

The customer is aware of being legally liable to Heidolph Instruments for any damages arising from incomplete and incorrect information.

Date _____ **Signature** _____

Company stamp

Please note

The shipper is responsible for the return of the goods in well packed condition, suitable for the mode of transport.

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This is a translation of the original operating manual. We reserve the right to make any technical modifications without prior notification. This documentation is not subject to revision; the latest version of this documentation can be found on our homepage. The copyright for all texts and images is held by Heidolph Instruments GmbH & Co. KG.
© Heidolph Instruments GmbH & Co. KG