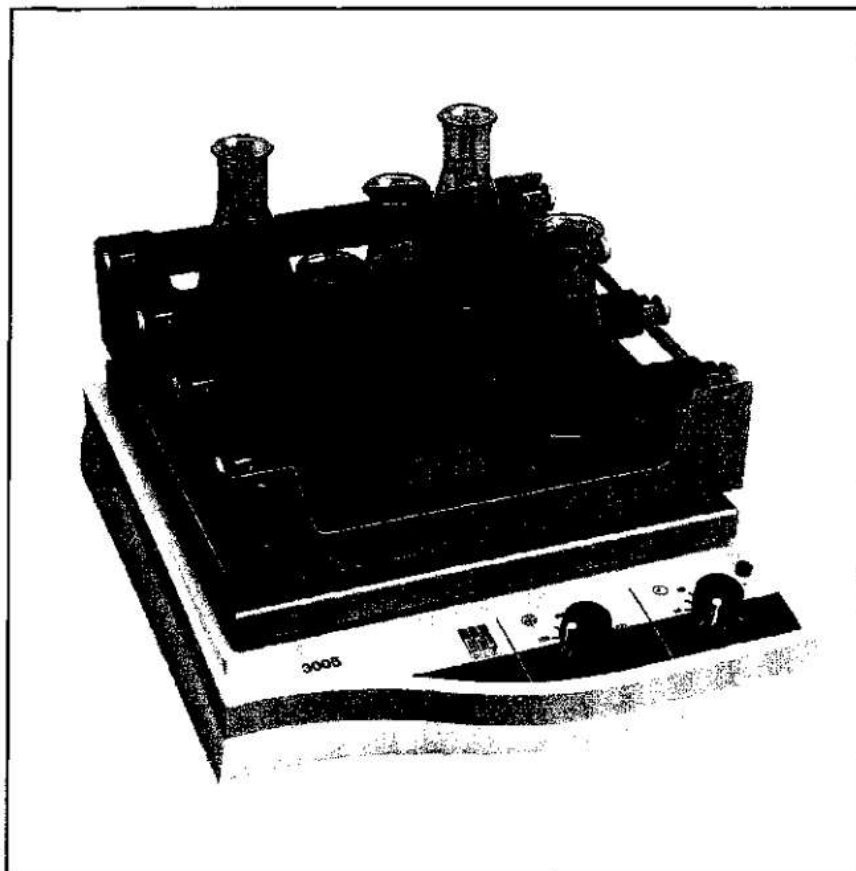
**Schüttelapparate
Shakers**

**3005-3006-3011-3013
3015-3016-3019-3023**

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**Operating
Instructions**

Schüttelapparate - Shakers

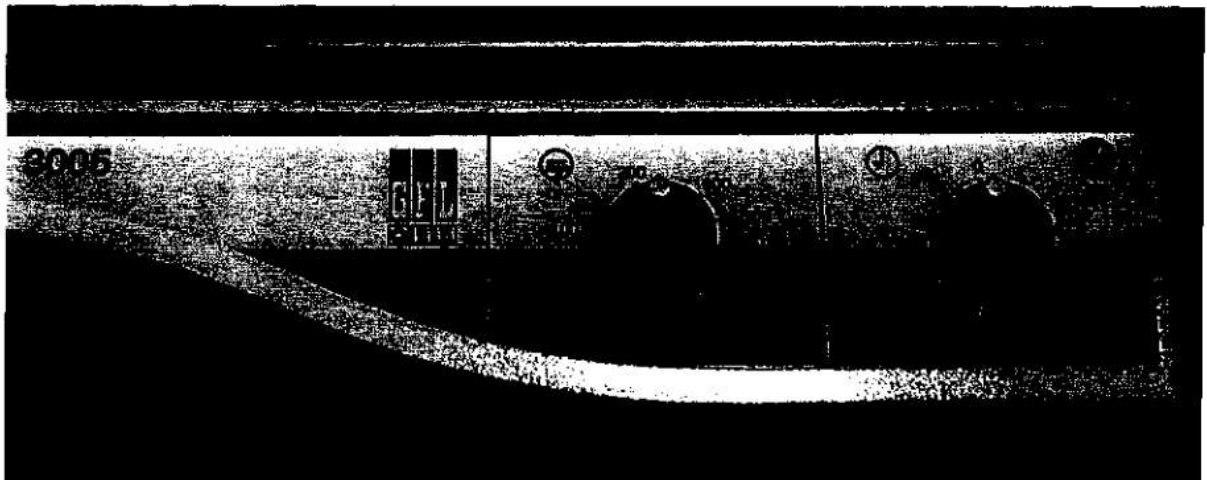
3005 / 3006 / 3011 / 3013 / 3015 / 3016 / 3019 / 3023

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Electronically geared AC motors as well as stable and durable shaking mechanisms create eight different Shaker types for four different motions: orbital, reciprocating, orbital rocking and rocking.

GFL Shakers are extremely silent and universally applicable, suitable for both gentle moving of liquids and vigorous mixing.

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Control panel

-to choose the operation
mode: continuous or timed
run

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Before installation, please check whether contents of package are in good order and complete.

Should you note any damages or have any reasons for complaint, please contact your supplier or directly:

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1. Before installation

The information in these operating instructions must **by all means** be carefully read and observed. Only then can a perfect operation of the **Shaker** be guaranteed.

Vital information within the manual are emphasized in bold letters. Safety precautions are additionally marked with the following symbols

Warning of dangerous electrical voltage

Warning of hand injuries

General hazard warning

A free of charge guarantee repair cannot be granted for defects due to improper installation or handling.

In some Shaker types, a paperboard strip has been inserted as a transport protection between shaking table and housing. The paperboard strip must be removed before installation.

2. Location of the Shaker

The unit is suitable for bench mounting **indoors** and for use in incubation and tempering rooms. Place on solid, stable and level surfaces only. Please observe the total admissible weight (weight of the unit as per Technical Data – item 7 of these operating instructions – plus load).

For use in incubation and tempering rooms please observe the environmental conditions stated in the Technical Data.

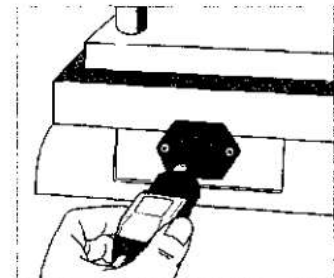
The Shaker is **not** suitable for use in **explosion endangered surroundings**, eg during anaesthesia with inflammable gas or steam types!

3. Voltage

The right-hand control knob of the **Shaker** must be in **0 position**.

The voltage on the nameplate (at the back of the unit) must be identical to the mains voltage.

If they are identical, connect the unit to the mains: Plug the appliance plug into the unit socket at the back of the Shaker and the shock-proof plug into the mains socket.

Schüttelapparate - Shakers

3005 / 3006 / 3011 / 3013 / 3015 / 3016 / 3019 / 3023

The shaker must only be connected to a properly installed power connection with neuter (PE), according to the local regulations. It must be secured that power can be cut all-pole (switch or socket).

4. Initiation

Caution,
increased danger of injuries during operation!
Do not reach into the unit during operation!

Caution,
by all means make sure to prevent the shaking procedure
creating an explosive atmosphere
during operation of the shaker!

Before initiation turn the left-hand control knob anti-clockwise to limit stop (minimum shaking frequency). Then turn right-hand control knob clockwise until the green pilot lamp glows.

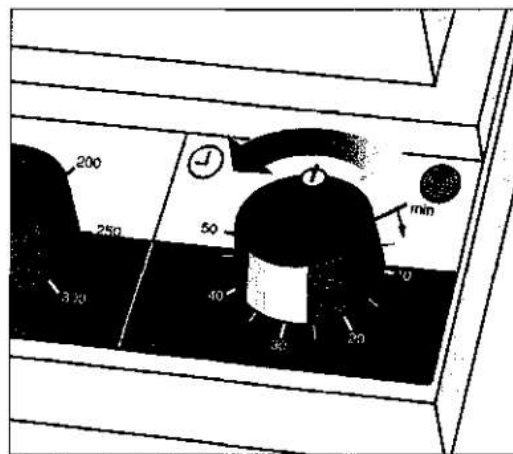
4.1 Choosing the operation mode

For **continuous operation** turn the knob to the first scale line; the knob will catch there. To switch off, turn the knob anti-clockwise to position 0.

For **timed runs** turn the control knob past the scale line for continuous run to the scale graduation marks for timed runs, showing time limits in minutes. The maximum timed period is 60 minutes. After expiry of the timed run, the Shaker will be switched off automatically.

Timed runs chosen by mistake can be reversed by turning the control knob anti-clockwise.



To **switch off** the Shaker, turn the control anti-clockwise to position 0.

4.2 Choosing the shaking frequency

Turn the left-hand control knob clockwise to increase the shaking frequency and anti-clockwise to decrease the shaking frequency.

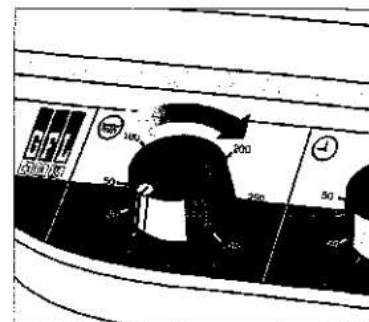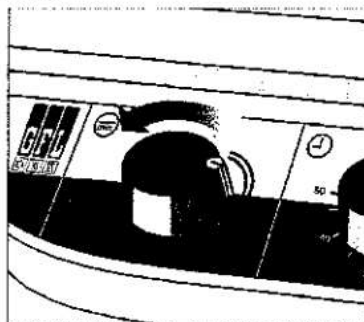The shaking frequency ranges of all Shaker types are stated in the Technical Data – item 7 – of these operating instructions.

5. Operation

A mechanical clockwork serves to run the Shaker in continuous or timed operation.

All models are equipped with ac motors, protected against overload. The electronic frequency control ensures a gentle start-up, irrespective of load.

A stable and durable mechanical construction serves to impart the torque of the motor with the specific shaking motion to the shaking platform.

6. Servicing, maintenance and clearing possible defects

Please make sure that no fluids come into contact with cable connections or the electrical parts of the inside unit!

Before cleaning and/or opening the unit (only by an electrician!) always pull the plug from the mains socket.

Danger of electrical shock !

Exchanging the fuses

The two fuses of the shaker are situated in a drawer below the unit plug. In order to check or to exchange these fuses, press the fixing clips on both sides of the drawer and pull.

The fuses may only be exchanged against fuses of the same type. Information on the fuse type can be found on the nameplate next to the unit plug and in the spare parts list of these operating instructions.

Breakdown in case of overload and mains failure

Overheating of the shaker due to overload will cause the shaker to be switched off.

Caution, the unit will restart automatically after cooling down or after mains failures.

If the unit is switched off due to any kind of breakdown, it must by all means be switched off, before the shaker is touched.

GFL Shakers are produced with first class materials and are made to withstand even rough service conditions. Nevertheless, the units should only be subjected to rough conditions within sensible limits.

The off-white pulver-coated surfaces of the housing and the shaking platform may be cleaned with mild detergents, if necessary.

Please do not hesitate to contact your local dealer or us for technical support.

In case of complaints please contact your local dealer or us:

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Servicings, repairs or modifications must be carried out according to the commonly recognised Technical Rules and Regulations by competent electricians only.

Only original spare parts must be used. Always demand a detailed confirmation of the carried out tasks by the person in charge (company, date, signature).

Schüttelapparate - Shakers

3005 / 3006 / 3011 / 3013 / 3015 / 3016 / 3019 / 3023

7. Technical data

Exterior dims. (wxdxh)	Model 3005	380 x 510 x 140mm / 330 x 330mm
Shaking platform (wxd)	Model 3006	380 x 510 x 140mm / 330 x 330mm
	Model 3011	510 x 625 x 170mm / 450 x 450mm
	Model 3013	510 x 625 x 170mm / 450 x 450mm
	Model 3015	510 x 625 x 142mm / 450 x 450mm
	Model 3016	510 x 625 x 145mm / 450 x 450mm
	Model 3019	745 x 730 x 135mm*/ 676 x 540mm
	Model 3023	380 x 510 x 125mm / 330 x 330mm
		* dimension without support rod
Max. load	Models 3005 and 3006	to max. 8 kg
> depending on the kind of load	Models 3011, 3013, 3015, 3016	to max. 15 kg
(even load distribution) and the	Model 3019	to max. 30 kg
required shaking frequency	Model 3023	to max. 1.2 kg
Motion /	Model 3005	orbital / 10mm / 20-500min ⁻¹
Motion amplitude /	Model 3006	reciprocating / 20mm / 20-300min ⁻¹
Motion frequency	Model 3011	three-dimensional orbital rocking / 3 degrees to horizontal / 2 – 50min ⁻¹
	Model 3013	rocking / 3 degrees to horizontal / 2 – 50min ⁻¹
	Model 3015	orbital / 30mm / 20-300min ⁻¹
	Model 3016	reciprocating / 30mm / 20-300min ⁻¹
	Model 3019	orbital / 32mm / 20-250min ⁻¹
	Model 3023	orbital / 3,0mm / 100-1450min ⁻¹
Electrical connection	all models	115V or 230V, +/-10%, 50....60Hz
Mains fuse	all models, house mains	circuit protected by 10 Amp
Power	Model 3005	82W
	Model 3006	40W
	Models 3011 and 3013	90W
	Models 3015, 3016 and 3023	65W
	Model 3019	90W
Protection type/class	all models	I / IP20
Surrounding conditions	all models	Use only inside buildings (not in explosion endangered surroundings)
Temperature	all models	+10°C to +50°C
Humidity	all models	max. 70% rel. humidity up to 31°C, decreasing to 50% rel. humidity at 50°C, non-condensing.
Weight	Models 3005 and 3006	9.1kg
	Model 3011	17.4kg
	Model 3013	15.4kg
	Models 3015 and 3016	16.8kg
	Model 3019	32.7kg
	Model 3023	12.5kg

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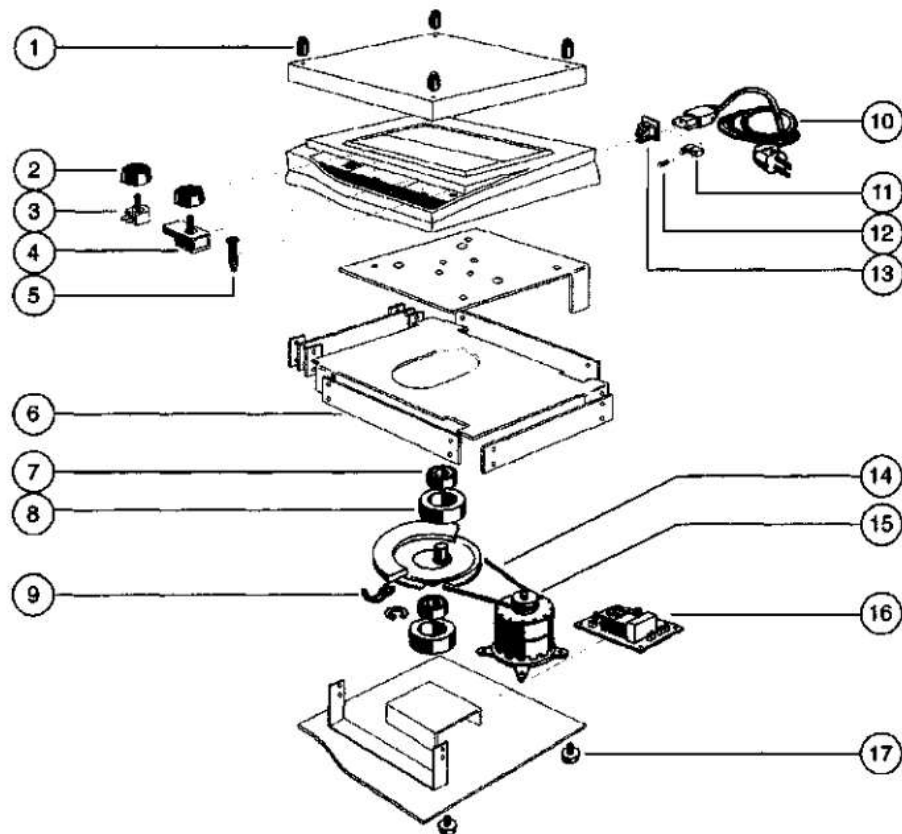
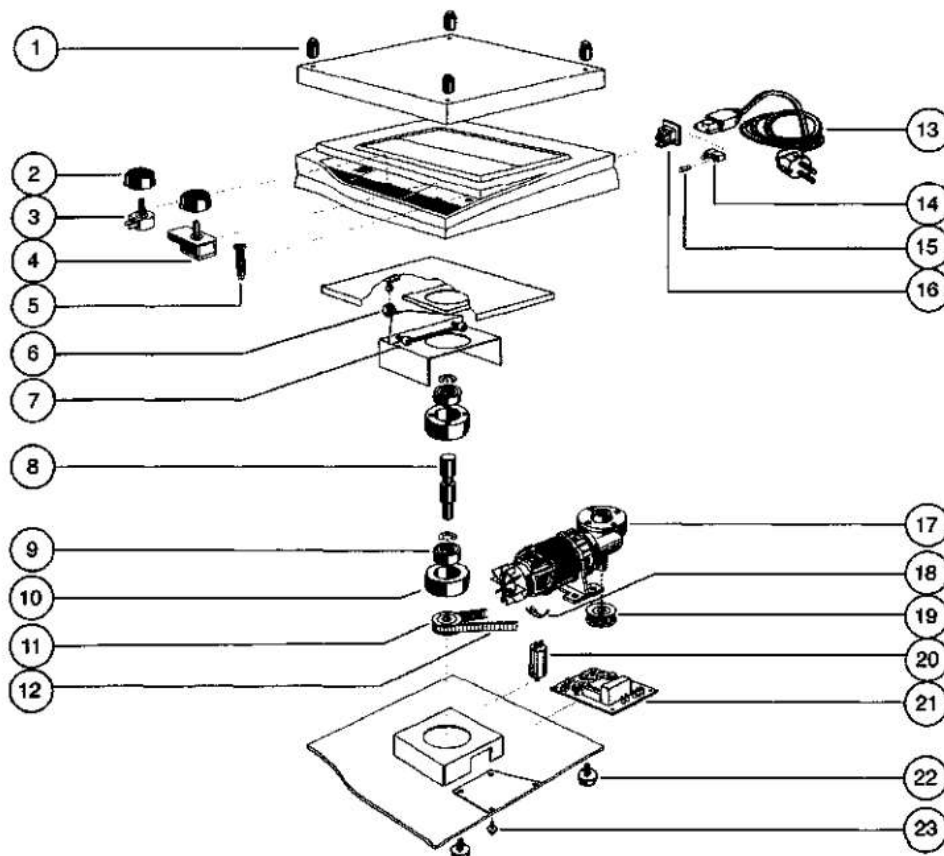
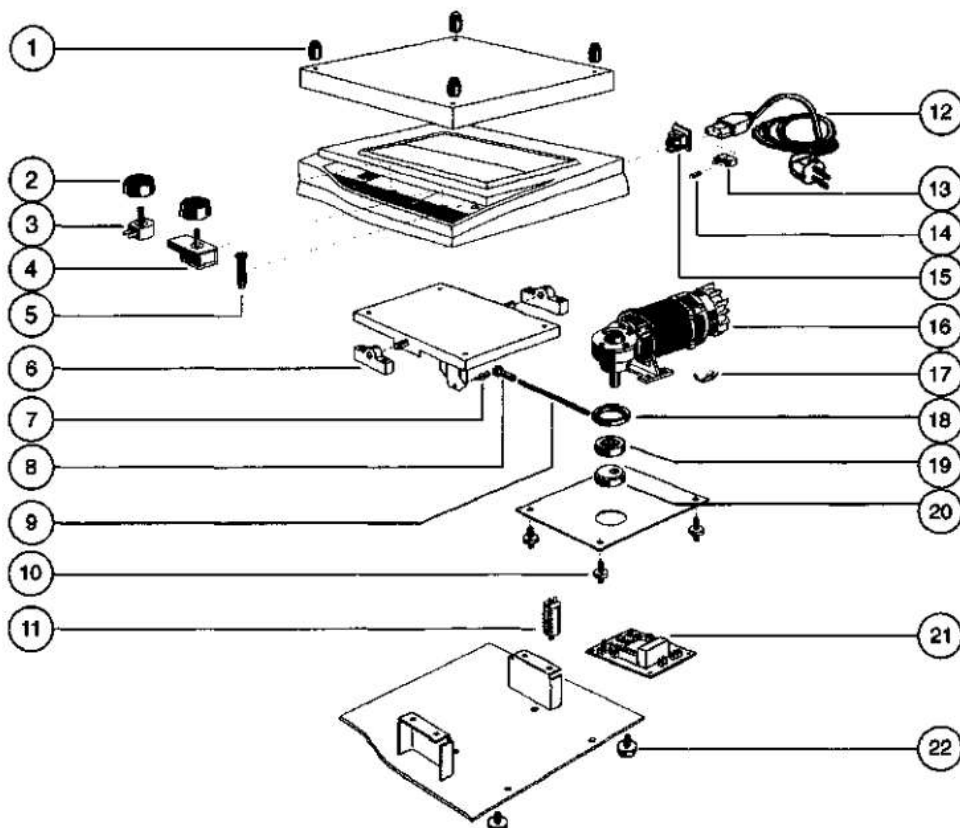
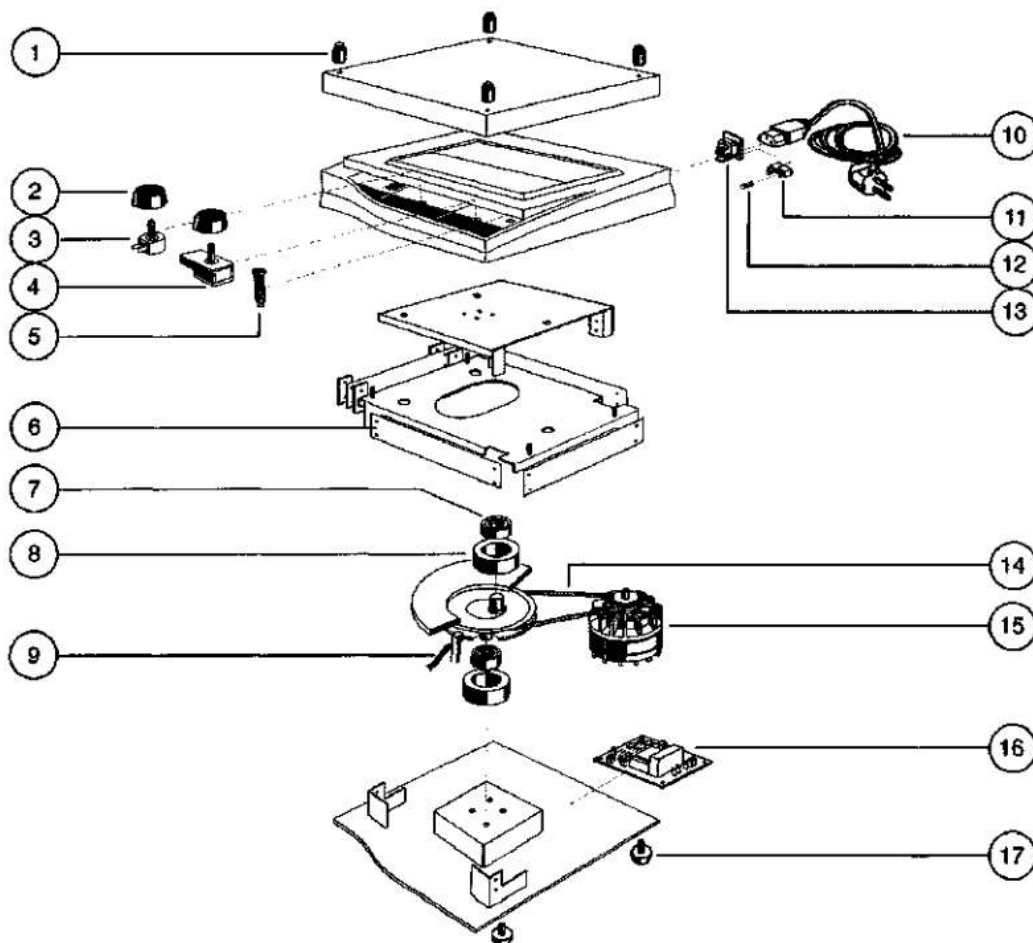
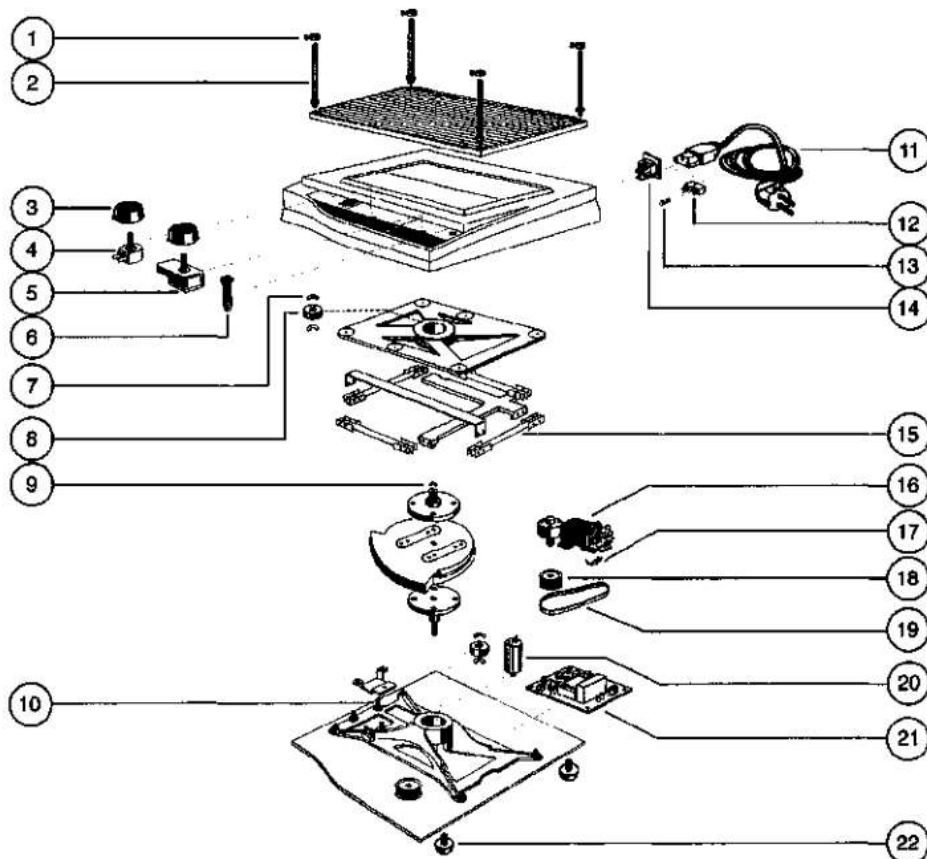
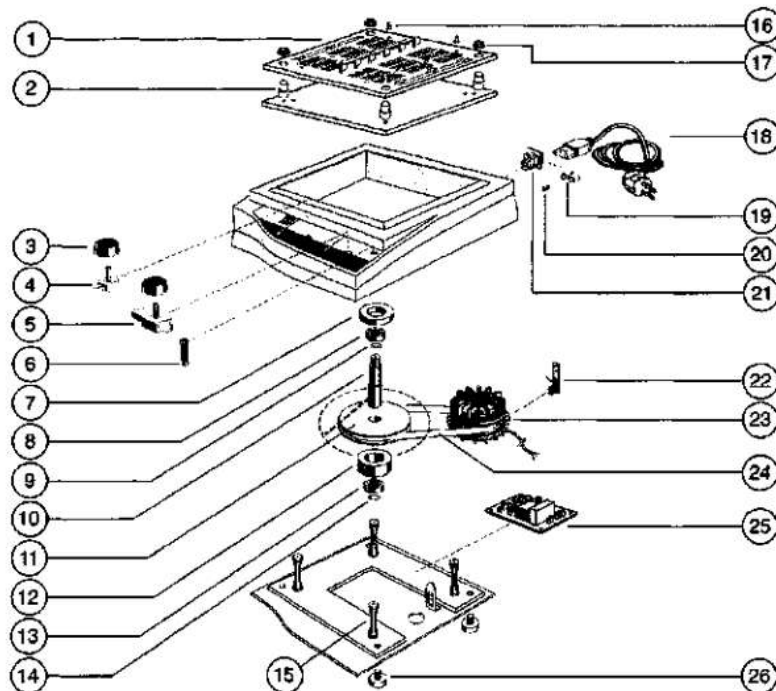
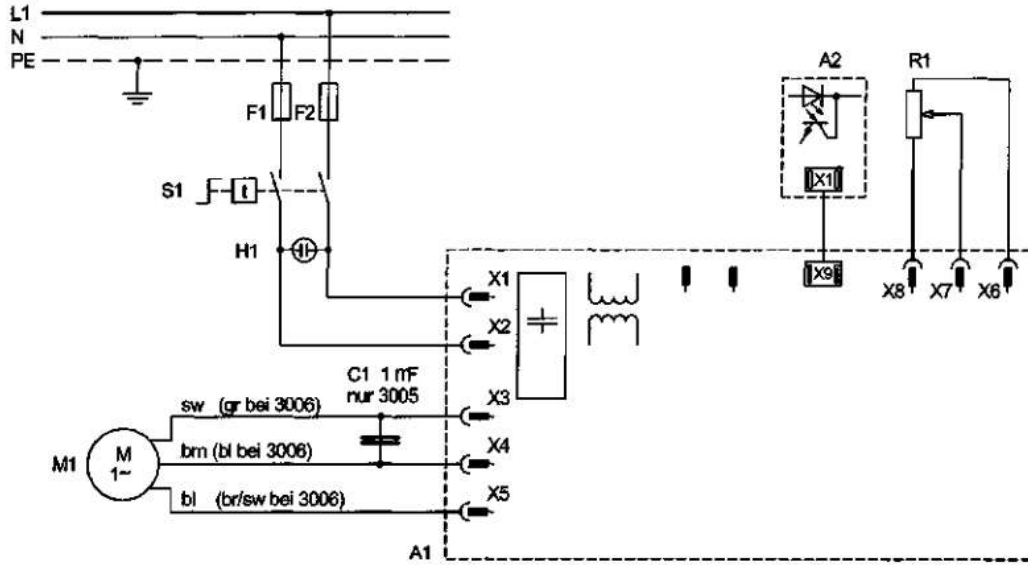
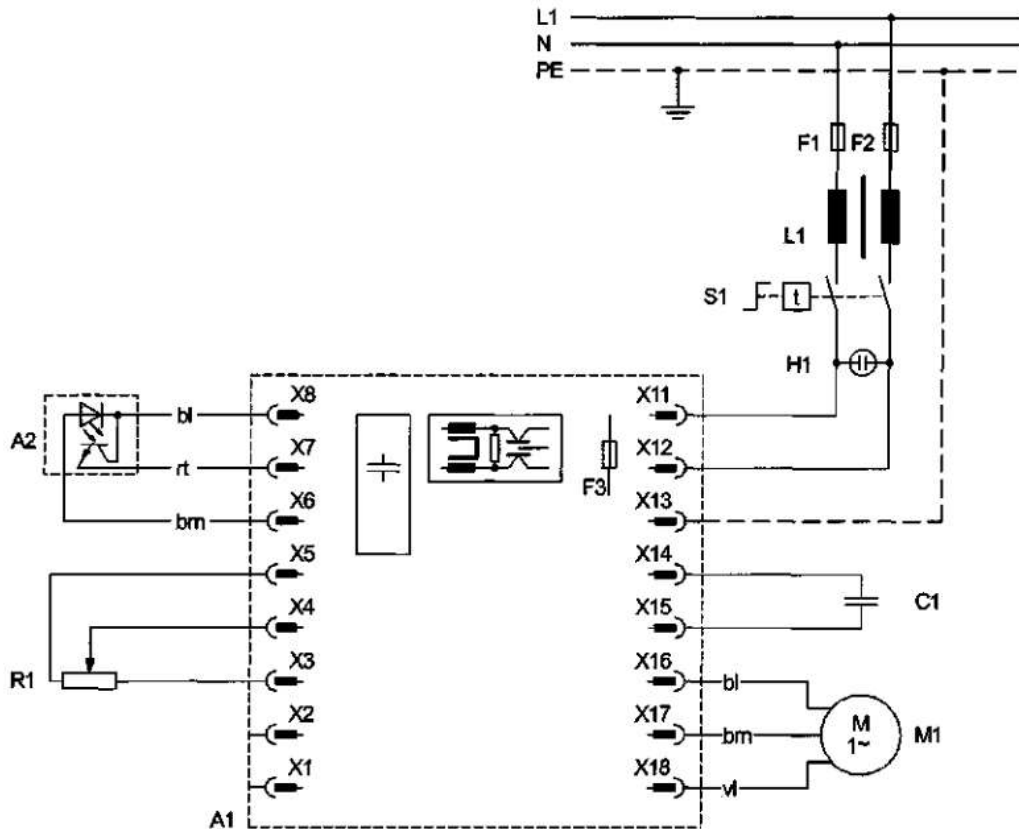
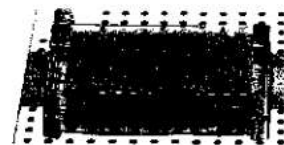
Order no. 3910
one set, including screwing
material

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Shaking Tray with holes to fix clamps for Erlenmeyer flasks and separating funnels, and to fix test tube rack 3953.

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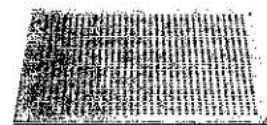
Order no. 3950,
made of stainless steel,
for models 3005 and 3006

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Order no. 3966,
made of stainless steel,
for models 3001, 3013, 3015
and 3016

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Order no. 3980,
made of anodized aluminium,
for model 3019

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Test Tube Rack, seat for test tubes can be tilted by 90° for easy loading, with springs for firm hold and silent shaking of the tubes. Supplied complete with fixing material to be screwed onto a shaking tray.

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Clamps for Erlenmeyer flasks, made of stainless steel. Supplied complete with fixing material to be screwed onto a shaking tray.

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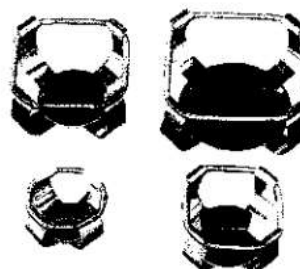
Clamps for separating funnels, made of stainless steel. Supplied complete with fixing material, to be screwed onto a shaking tray.

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Universal mount for safe fixing of various shaking objects between rubber-coated bars.

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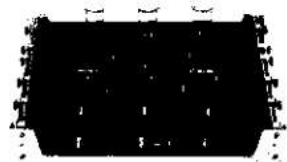
Platform Frames to increase the capacity.

Bestell-Nr.3968

für Typen 3011 und 3013

aus "Edelstahl Rostfrei"

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Oder no. 3968

for models 3011 and 3013,

made of stainless steel,

with 4 platforms.

Bestell-Nr.3981

für Typ 3019

aus "Edelstahl Rostfrei" zur

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3980

Order no. 3981

for model 3019,

made of stainless steel, to

accept 2 trays 3980.